



# VARIAN

## Application Note SI-01622

### SEC Analysis of Xanthan

Greg Saunders  
Varian, Inc.

#### Introduction

Xanthans are high molecular weight polymers that have the ability to raise the viscosity of water dramatically. This is exploited in oil recovery processes, where xanthan acts as a water thickener. Analysis by SEC is possible at very low sample concentrations, however, to produce significant detector response, an increased injection volume is required. The sample was analyzed using PL aquagel-OH 60 15  $\mu\text{m}$  columns, which were employed in order to avoid on-column shear degradation of these high molecular weight samples, and cover a molecular weight range from  $10^5$  to  $10^7$ .

#### Conditions

Sample: Xanthan  
Columns: 2 x PL aquagel-OH 60 15  $\mu\text{m}$ , 300 x 7.5 mm  
Eluent: 0.2 M  $\text{NaNO}_3$  + 0.01 M  $\text{NaH}_2\text{PO}_4$  at pH 7  
Flow Rate: 1.0 mL/min  
Detection: RI

#### Conclusion

SEC using PL aquagel-OH 60 15  $\mu\text{m}$  columns successfully analyzed a sample of xanthan at a very low concentration. Aqueous SEC with PL aquagel-OH columns provides information not only on the molecular weight of the polymer but also on the polydispersity and the shape of the molecular weight distribution. The excellent chemical and mechanical stability of these columns offer high performance with good repeatability and column lifetime.

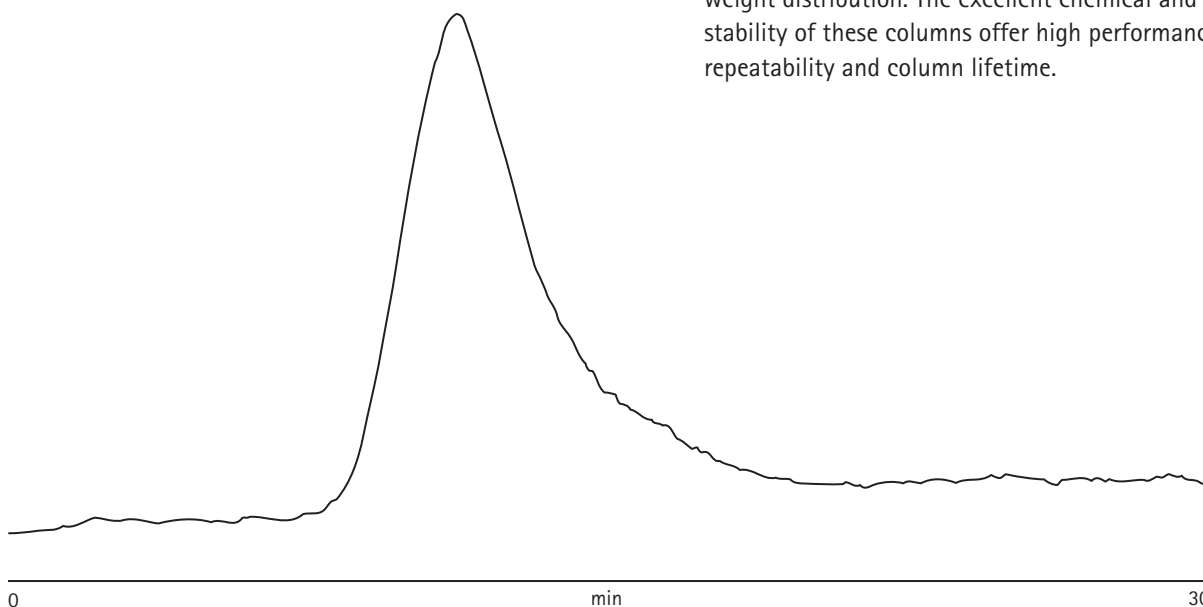


Figure 1. Raw data chromatogram of xanthan.

*These data represent typical results. For further information, contact your local Varian Sales Office.*

*Varian and the Varian Logo are trademarks or registered trademarks of Varian, Inc. in the U.S. and other countries.  
© 2009 Varian, Inc.*

Varian, Inc.  
www.varianinc.com  
North America: 800.926.3000 – 925.939.2400  
Europe: *The Netherlands*: 31.118.67.1000  
Asia Pacific: *Australia*: 613.9560.7133  
Latin America: *Brazil*: 55.11.3238.0400



# VARIAN